



## Quick Start Guide

---

Kima-Connect: Tubing Set for Kima™ Pump

## Kima-Connect: Tubing Set for Kima Pump

Kima-connect: tubing set for the Kima pump contains:

- 2 pump priming cables
- 1 pump inlet cable
- 1 pump outlet cable
- 1 biochip outlet cable

To connect the tubing to the Kima pump, follow the steps illustrated below.




Remove cover from pump. Connect pump priming cables to inlet and outlet of pump and prime with solution (e.g. deionised water or PBS) using the syringe.



Connect inlet cable as shown



Replace cover on pump. Connect pump outlet cable as shown.

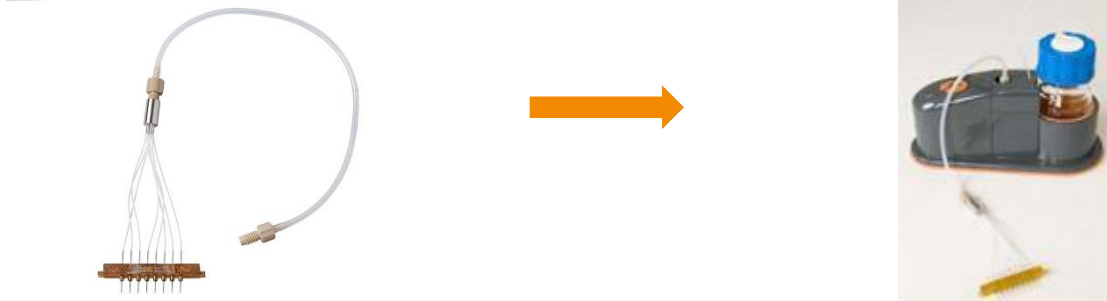

Connect biochip outlet cable as shown.

